illumina

Infinium[®] HumanCore-24 v1.2 BeadChip

Customizable array for cost-effective large-scale genotyping and screening studies.

Overview

Customizable Infinium HumanCore-24 v1.2 BeadChips offer an economical way to perform and support large genetic studies, especially large-scale genotyping studies. Developed in collaboration with several leading research institutions, HumanCore-24 BeadChips contain more than 240,000 highly informative genome-wide tag single nucleotide polymorphisms (SNPs) and over 20,000 high-value markers, including insertions/deletions (indels) and updated exomefocused content (Table 5 and Table 6). The Infinium HumanCore-24+ v1.2 BeadChip has the added capacity to include up to 300,000 semicustom markers. In addition to performing cost-effective largescale genotyping studies, HumanCore-24 BeadChips can be used to obtain baseline sample data sets for various downstream applications quickly and easily. These applications include common variant, mitochondrial DNA (mtDNA), ancestry, sex confirmation, loss ofvariant, and indel, detection studies. Infinium HumanCore-24 v1.2 BeadChips are based on the trusted Infinium Assay. Using the proven iScan[™] or HiScan[™] System, integrated analysis software, and Infinium high-throughput screening (HTS) Assay, this high-density, 24-sample BeadChip (Figure 1) combines affordability with high-throughput sample processing to deliver high-quality, genome-wide information.

			_
<u> </u>	_		_
	_		
			-
		_	_
			_
	=		_
	_		
		_	
			_
		_	
	illun	nina	
02	7416002	20.24-1	GP
93	7418003	J7 Z4X -	AD.
100			III III

Figure 1: The Infinium HumanCore-24 v1.2 BeadChip— The Infinium HumanCore-24 v1.2 BeadChip enables genotyping of markers informative across diverse populations, delivering high-quality data that can be used for various downstream applications.

High-throughput workflow

The Infinium HumanCore-24 v1.2 BeadChip uses the highly scalable 24-sample Infinium HTS format for high-throughput processing of thousands of samples per week for large, population-scale research and variant screening. The Infinium HTS format also provides a rapid three-day workflow that allows genotyping service providers and clinical researchers to gather data and advance studies quickly (Figure 2).

Optional integration of the Illumina Laboratory Information Management System (LIMS) into the workflow provides high laboratory efficiency with automation functionality, process tracking, and quality control (QC) data tracking. The Illumina ArrayLab Consulting Service offers customized solutions to high-throughput genotyping labs that desire increased efficiency and overall operational excellence.

Robust, high-quality assay

The Infinium HumanCore-24 v1.2 BeadChip uses proven Infinium assay chemistry to deliver the same high-quality, reproducible data (Table 1) that Illumina genotyping arrays have provided for over a decade. The Infinium product line provides high call rates and high reproducibility for numerous sample types including, saliva, blood, solid tumors, fresh frozen, and buccal swabs (Table 2–4). In addition, the high signal-to-noise ratio of the individual genotyping calls from the Infinium assay provides researchers with access to genome-wide copy number variant (CNV) calling with a mean probe spacing of ~ 9.5 kb.

Design Markers | Process BeadChip | Analyze Data



Figure 2: The Infinium HTS workflow–The Infinium HTS format provides rapid 3day workflow with minimal hands-on time.

Table 1: Product information

Table 1. Floudet mornation				
Feature	Description			
Species	Human			
Total number of markers	305,445			
Capacity for custom bead types	300,000			
Number of samples per BeadChip	24 Samples			
DNA input requirement	200 ng			
Assay chemistry	Infinium HTS			
Instrument support	iScan or HiScan System			
Sample throughput ^a	~ 2304 samples/week			
Coop time ner comple	iScan System	HiScan System	ו	
Scantime per sample	2.5 min	2.0 min		
Data performance	Value ^b	Product Spec	ificationd	
Call rate	99.70%	> 99% avg.		
Reproducibility	99.99%	> 99.9%		
Log R deviation	0.09	< 0.30°		
Spacing				
Specing (kb)	Mean	Median	90th%°	
opacitily (ku)	9.5	5.8	21.7	
a Estimate essumes 1 iCoop I	Queteres 1 Autol	ander O.V. O.Te	aan valaata and	

 Estimate assumes 1 iScan System, 1 AutoLoader 2.x, 2 Tecan robots, and a 5-day work week.

b. Values are derived from genotyping 333 HapMap reference samples.

c. Value expected for typical projects using standard Illumina protocols. Tumor samples and samples prepared by methods other than standard Illumina protocols are excluded.

d. Excludes Y chromosome markers for female samples.

Table 2: Imputation accuracy from 1000G^a at various MAF thresholds

Deputation ^b	I	mputation Accurac	y
Population	$MAF \ge 5\%$	$MAF \ge 1\%$	MAF 1-5%
AFR	0.90	0.84	0.76
AMR	0.94	0.89	0.79
EAS	0.92	0.85	0.66
EUR	0.94	0.89	0.76
SAS	0.92	0.86	0.70

a. Compared against Phase 3, version 5 of the 1000 Genomes Project (1000G). www.1000genomes.org. Accessed July 2016.

b. See www.1000genomes.org/category/frequently-asked-questions/population

Abbreviations: MAF: minor allele frequency; AFR: African; AMR: Ad Mixed American; EAS: East Asian; EUR: European; SAS: South Asian.

Table 3: LD $r^2 \ge 0.80$ from 1000G^a at various MAF thresholds

1000C Dopulation ^b	LD Coverage ($r^2 \ge 0.80$)		
rooud Population	$MAF \ge 5\%$	$MAF \ge 1\%$	
AFR	0.27	0.16	
AMR	0.55	0.38	
EAS	0.65	0.53	
EUR	0.61	0.47	
SAS	0.57	0.42	

a. Compared against Phase 3, version 5 of the 1000 Genomes Project (1000G). www.1000genomes.org. Accessed July 2016.

b. See www.1000genomes.org/category/frequently-asked-questions/population

Abbreviations: LD: linkage disequilibrium; AFR: African; AMR: Ad Mixed American; EAS: East Asian; EUR: European; SAS: South Asian.

Table 4: LD mean r² from 1000G^a at various MAF thresholds

Dopulation ^b	LD Coverage (Mean r ²)		
Fopulation	MAF ≥ 5%	$MAF \ge 1\%$	
AFR	0.45	0.28	
AMR	0.70	0.50	
EAS	0.76	0.62	
EUR	0.73	0.57	
SAS	0.71	0.54	

a. Compared against Phase 3, version 5 of the 1000 Genomes Project (1000G). www.1000genomes.org. Accessed July 2016.

b. See www.1000genomes.org/category/frequently-asked-questions/population

Abbreviations: LD: linkage disequilibrium; MAF: minor allele frequency; AFR: African; AMR: Ad Mixed American; EAS: East Asian; EUR: European; SAS: South Asian.

Table 5: Marker information

Marker categories			No. of markers
Exonic markers ^a			41,698
Intronic markers ^a			120,454
Nonsense markers ^b			9833
Missense markers ^b			6707
Synonymous markers ^b			5430
Mitochondrial markers ^c			161
Indels ^c			12,312
Say abromage	Х	Y	PAR/homologous
Sex chromosomes	8100	2004	152

a. RefSeq - NCBI Reference Sequence Database.

www.ncbi.nlm.nih.gov/refseq. Accessed September 2016.b. Compared against the University of California, Santa Cruz (UCSC) Genome Browser. genome.ucsd.edu. Accessed August 2014.

c. NCBI Genome Reference Consortium, Version GRCh37.

www.ncbi.nlm.nih.gov/grc/human. Accessed July 2016.

Abbreviations: indel: insertion/deletion; PAR: pseudoautosomal region.

Table 6: High-value content

Content	No. of markers	Research application/note
ADME core and extended genes ¹	5904	Drug metabolism and excretion
Blood phenotype genes ³	255	Blood phenotypes
COSMIC ⁴ genes	137,811	Somatic mutations in cancer
GO ⁵ CVS genes	37,104	Cardiovascular conditions
Database of genomic variants ⁶	235,945	Genomic structural variation
eQTLs ⁷	2438	Genomic loci regulating mRNA expression levels
Fingerprint SNPs ⁸	127	Human identification
HLA Genes ²	60	Disease defense, transplant rejection, and autoimmune disorders
Extended MHC ^{a9}	1430	Disease defense, transplant rejection, and autoimmune disorders
KIR genes ²	7	Autoimmune disorders and disease defense
Neanderthal SNPs ¹⁰	352	Neanderthal ancestry and human population migration
NHGRI GWAS Catalog ¹¹	5158	Markers from published genome-wide association studies
RefSeq ¹² 3' UTRs	8541	3' untranslated regions of known genes
RefSeq 5' UTRs	3691	5' untranslated regions of known genes
RefSeq All UTRs	11,890	All untranslated regions of known genes
RefSeq	154,358	All known genes
RefSeq +/- 10 kb	177,866	All known genes plus regulatory regions
RefSeq Promoters	6345	2 kb upstream of all known genes to include promoter regions
RefSeq splice regions	5325	Variants at splice sites in all known genes

a. Extended MHC is a ~ 8 Mb region.

Abbreviations: ADME: absorption, distribution, metabolism, and excretion; APOE: apolipoprotein E; COSMIC: catalog of somatic mutations in cancer; GO CVS: gene ontology annotation of the cardiovascular system; eQTL: expression quantitative trait loci; HLA: human leukocyte antigen; KIR: killer cell immunoglobulin-like receptor; MHC: major histocompatibility complex; NHGRI: national human genome research institute; GWAS: genome-wide association study; UTR: untranslated region; RefSeq: reference sequence.

Ordering information

Infinium HumanCore-24 v1.2 Kit	Catalog no.		
48 samples	20024566		
288 samples	20024567		
1152 samples	20024568		
Infinium HumanCore-24+ v1.2 Kit ^a	Catalog no.		
48 samples	20024569		
288 samples	20024660		
1152 samples	20024661		
a. Enabled for additional custom content.			

Learn more

To learn more about the Infinium HumanCore-24 v1.2 BeadChip and other Illumina genotyping products and services, visit www.illumina.com/genotyping

References

- 1. PharmaADME Gene List. www.pharmaadme.org. Accessed August 2014.
- 2. University of California, Santa Cruz (UCSC) Genome Browser. genome.ucsc.edu. Accessed August 2014.
- NCBI Reference Sequence Blood Group Antigen Gene Mutation Database. www.ncbi.nlm.nih.gov/projects/gv/rbc/xslcgi.fcgi?cmd=bgmut/systems. Accessed July 2016.
- 4. Catalog of somatic mutations in cancer. cancer.sanger.uk/cosmic. Accessed July 2016.
- 5. Gene Ontology Consortium. www.geneontology.org. Accessed July 2016.
- Database of Genomic Variants. dgv.tcag.ca/dgv/app/home. Accessed July 2016.
- NCBI eQTL Database. www.ncbi.nlm.nih.gov/projects/gap/eqtl/index.cgi. Accessed July 2016.
- 8. The Allele Frequency Database. alfred.med.yale.edu/alfred/snpSets.asp. Accessed July 2016.
- de Bakker PIW, McVean G, Sabeti PC, et al. A high-resolution HLA and SNP haplotype map for disease association studies in the extended human MHC. *Nat Genet*. 2006;38:1166–1172.
- Neanderthal Genome Browser. neandertal.ensemblgenomes.org/index.html. Accessed July 2016.
- 11. NHGRI GWAS Catalog. www.ebi.ac.uk/gwas/docs/downloads. Accessed July 2016.
- 12. NCBI Reference Sequence Database. www.ncbi.nlm.nih.gov/refseq. Accessed July 2016.